
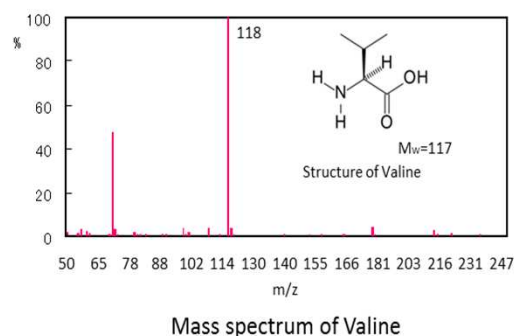


## アミノ酸の分離 (LC/MS)

### Amino acids (LC/MS)

Sunniest RP-AQUA 5  $\mu$ m, 150 x 2.1 mm i.d.

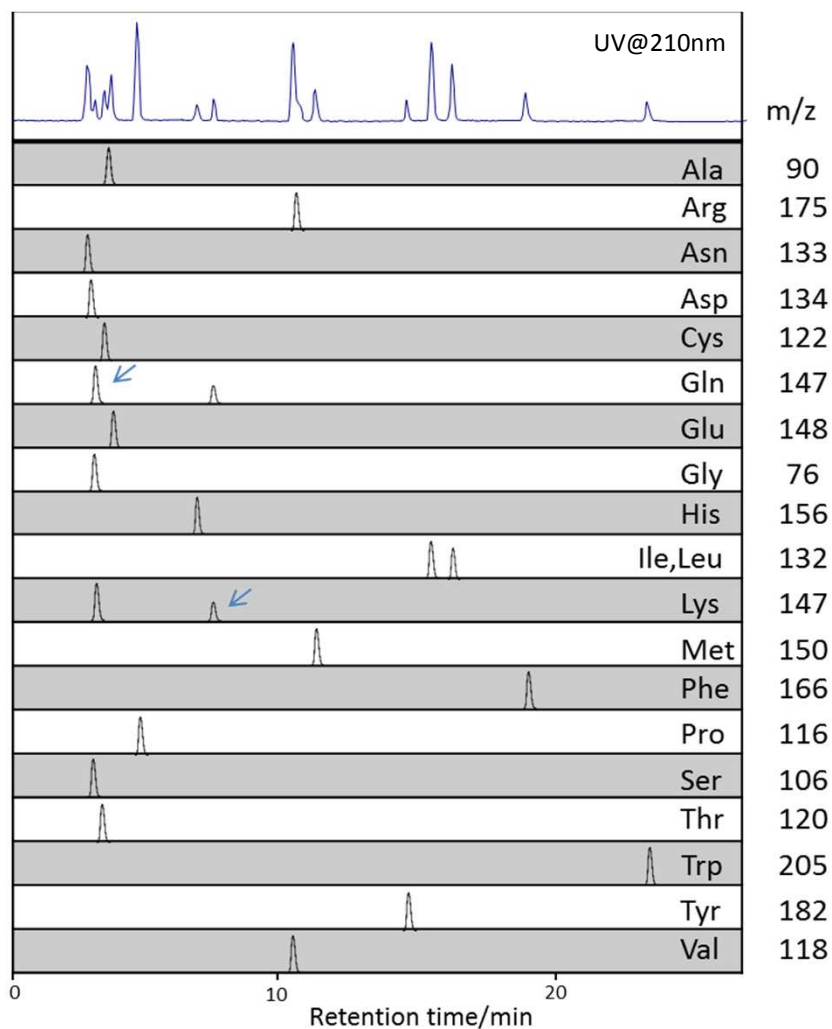

Column: Sunniest RP-AQUA 5  $\mu$ m, 150 x 2.0 mm

Mobile phase:

A) 5mM HFBA (Heptafluorobutyric acid)

B) 5mM HFBA in Acetonitrile/water(9/1)

%B 0% to 20% in 20 min

Flow rate: 0.2mL/min

Temperature: 40  $^{\circ}$ C

Detection: Quattro Micro API (ESI positive) SIM and UV@210nm