
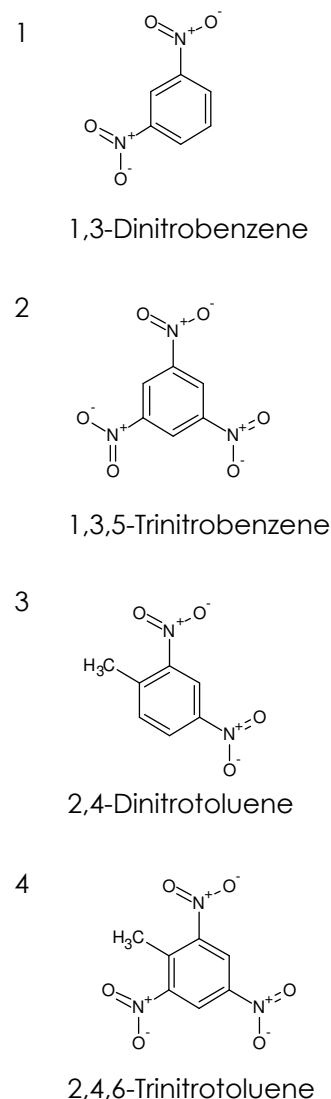


爆薬成分の分離 (2)

Components of explosive(2)

SunShell C18 2.6 μm, 100 x 4.6 mm i.d.
 SunShell PFP 2.6 μm, 100 x 4.6 mm i.d.
 SunShell RP-AQUA 2.6 μm, 100 x 4.6 mm i.d.
 SunShell Phenyl 2.6 μm, 100 x 4.6 mm i.d.
 SunShell Biphenyl 2.6 μm, 100 x 4.6 mm i.d.

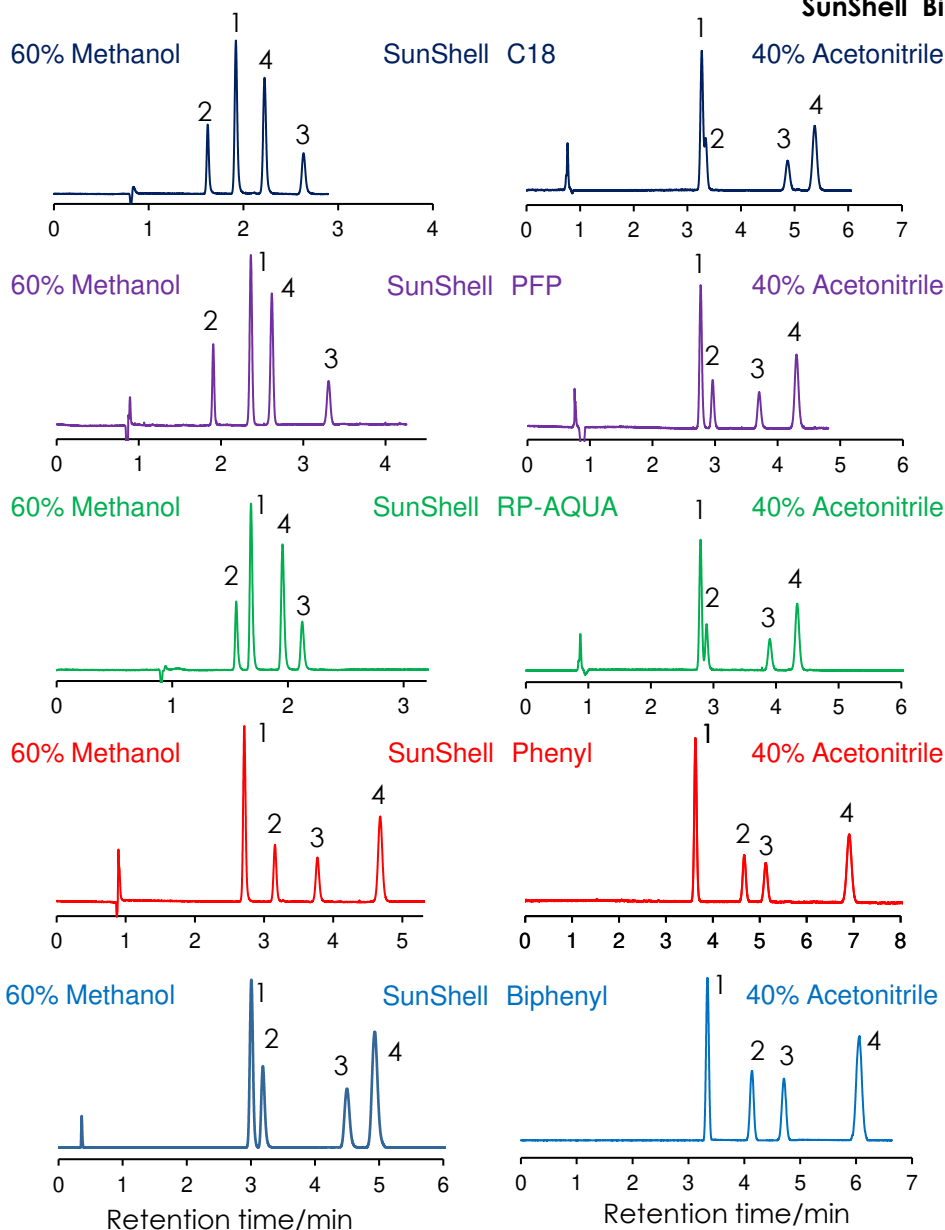

Column: SunShell C18 2.6 μm, 100 x 4.6 mm
 SunShell PFP 2.6 μm, 100 x 4.6 mm
 SunShell RP-AQUA 2.6 μm, 100 x 4.6 mm
 SunShell Phenyl 2.6 μm, 100 x 4.6 mm
 SunShell Biphenyl 2.6 μm, 100 x 4.6 mm

Mobile phase: Methanol:water=60:40
 Acetonitrile:water=40:60

Flow rate: 1.0 mL / min

Temperature: 40 °C

Detection: UV@250 nm

Sample: 1 = 1,3-Dinitrobenzene, 2 = 1,3,5-Trinitrobenzene, 3 = 2,4-Dinitrotoluene, 4 = 2,4,6-Trinitrotoluene